PRODUCT INFORMATION Colloidal Gold Conjugate

Catalog Number:	GP-2002-15
Description:	Colloidal Gold Conjugated <i>Ricinus communis</i> Lectin (Castor Bean) -RCA-II-, 15nm Particle Size, 1mL
Lot Number:	
Expiration Date:	
Protein Concentration:	
HauCl ₄ Concentration:	OD 520 = 4
Buffer:	Distilled H ₂ O, pH 7.0 - 7.5
Storage:	<u>DO NOT FREEZE</u> ! Store at 5-8°C.
Other Data:	Absorbance peak at 520nm. CONTAINS 0.02% SODIUM AZIDE AS A PRESERVATIVE. Gently resuspend any sediment. If necessary, clarify by centrifugation at 400-500 x g for ten minutes before use.
Quality Control:	Procedure developed by EY Laboratories, Inc.
References:	 Roth, J. Techniques in Immunocytochemistry. Eds. Bullock and Petrusz Academic Press 1983, 217-284.
CAUTION:	RCA-II is toxic. Decontaminate all glassware with bleach (5%).





Tel: 650-342-3296 Fax: 650-342-2648 Orders: 1-800-821-0044 (Outside CA only)

"TOXIC" MATERIAL SAFETY DATA SHEET

Effective Date: March 31, 2006 Revision 6 Page 1 of 2

PRODUCT IDENTIFICATION

Name: Ricin Alpha Chain and Ricin Beta Chain. Crude and purified Ricinus communis, Viscum album, Abrus precatorius. Also, these proteins conjugated to fluorescein isothiocvanate, tetramethylrhodamine isothiocvanate, Texas Red®, D-Biotin, Horseradish Peroxidase, Alkaline Phosphatase, and Ferritin. Catalog Number (s): AAL-2001, AL-2001, AL-2003, BA-2001, BA-2002, BA-4001, BA-5401, BAL-2001, F-2001, F-2002, F-4001, F-5401, FAL-2001, GP-2001, GP-2002, GP-4001, GP-5401, H-2001, H-2002, H-4001, H-5401, HAL-2001, I-2001, I-2002, I-4001, I-5401, L-2000, L-2001, L-2002, L-4000, L-4001, L-5400, L-5401, LA-2001, LA-2002, LA-4001, LA-5401, MB-2001, PB-2001, R-2001, R-2002, R-4001, R-5401, RAL-2001, T-2001, T-2002, T-4001, T-5401, TAL-2001, A-2001, A-2002, A-4001, A-5401, AK-2001, AK-2002, AK-4001, AK-5401, LGK-001, LGK-001A, LGK-002, LGK-002A, LGK-003, RAC-1, RBC-1. Formula: Complex polypeptides. Synonyms: RCA-I, RCA-II, RCA120, RCA60, castor bean lectin, ricin, ricin toxin, VAA, mistletoe lectin, viscumin, APA, abrin, jequirity bean lectin. These proteins may be

mistletoe lectin, viscumin, APA, abrin, jequirity bean lectin. These proteins may be coupled to various fluorescent dyes, enzymes, ferritin, or D-biotin. CAS Number: 1393-62-0 (for Abrus precatorius).

EMERGENCY INFORMATION

EY Laboratories, Inc. 107 North Amphlett Blvd. San Mateo, CA 94401

EFFECTS OF

ROUTES

EXPOSURE

OVEREXPOSURE

EMERGENCY PHONE: 650-342-3296

HAZARDOUS COMPONENTS

Specific protein (s) as listed on the vial label. These proteins are always supplied in solution, never in dried form. The concentration of the specific protein in solution generally will be greater than 0.1% (w/v, 1mg/1ml) and less than 1% (w/v, 10mg/ml) unless otherwise indicated on the vial label and the product information sheet. These proteins are toxins which function by inhibiting protein synthesis. Extreme caution should be used when handling these products. The fluorescent molecules fluorescein and tetramethylrhodamine isothiochyanate, as well as Texas Red® are possible carcinogens. Although the complete properties of these molecules have not been fully investigated, similar compounds are known to be chemically reactive with proteins and other biomolecules. Ferritim may be tumorigenic and should not be allowed to come in contact with skin. Alkaline Phosphatase, Horseradish Peroxidase, and D-Biotin may be harmful if ingested, inhaled, or allowed to contact the skin. Some individuals may exhibit a stronger adverse reaction to these chemicals than other individuals. All of the above toxins and related products should be handled only by qualified personnel and should be used in a facility isolated from the common work area. Sodium azide is included in most solutions as a preservative at a concentration of 0.02-1%.

HEALTH HAZARD INFORMATION

EXPOSURE LIMITS: No exposure limits have been established, however, no exposure is considered safe. Ricin (RCA-II, RCA60) has been tested for toxicity. The LD50 (iv) in mice is 2µg/100 grams body weight. The other toxins are less potent but specific toxicity

> And contact with these toxins may result in nausea, cramps, fatigue, and other flulike symptoms. Chronic or high dose exposure to the toxins may result in serious illness or death. These toxins function by inhibiting protein synthesis. Any living tissue may be affected by these toxins.

> Contact with skin and mucous membranes are the primary routes of exposure. Care should be taken to avoid the formation of aerosols when handling any of the solutions. The use of glass equipment is discouraged since sharp edges may puncture protective equipment and directly inject toxic material.

LABORATORIES, INC.

07 North Amphlett Blvd. San Mateo, CA 94401

```
Tel: 650-342-3296
Fax: 650-342-2648
Orders: 1-800-821-0044
(Outside CA only)
```

TOXIC MSDS for Crude and Purified Proteins and Enzymes Continued - page 2 of 2.

PHYSICAL CHARACTERISTICS

 APPEARANCE:
 Liquids will be clear to dark purple.

 SOLUBILITY:
 All liquids are completely miscible in water and buffers.

FIRE AND EXPLOSION HAZARDS

 Not considered to be a fire hazard. High heat completely will destroy any toxic activity of the proteins.

 EXTINGUISHING MEDIA:
 Water spray or CO₂.

 SPECIAL FIRE FIGHTING PRECAUTIONS:
 As required for the surrounding fire.

NOTE: Most solutions contain 0.02-1% sodium azide as a preservative. Azide may react with lead and copper plumbing to form explosive metal azides. Flush with copious amounts of water when disposing material in the sink.

REACTIVITY DATA

STABILITY:	Stable. Decomposition products are not known to be hazardous.
HAZARDOUS POLYMERIZATION:	Will NOT occur.
INCOMPATIBILITY:	None known. (Lead and copper may react with sodium azide).

SPILL / LEAK PROCEDURES

MATERIAL RELEASE / SPILL: Avoid contact with liquid. Wear a self-contained breathing apparatus and a full protective suit when cleaning up large spills. Flush contaminated area with a 0.5-5% solution of hypochlorite (household bleach diluted 1:5 with water is sufficient. Clean up spill with a paper towel soaked in household bleach. Do not allow solutions to dry on environmental surfaces. Wash affected area with detergent after the area has been treated with bleach.
 WASTE DISPOSAL: Incinerate, autoclave, or dispose of paper waste in accordance with all Local, State, and Federal regulations. All laboratory equipment and supplies that contact these toxins must be soaked in bleach for at least 1 hour or autoclaved. All of these proteins are generally not considered environmental hazards once treated with bleach.

EMERGENCY FIRST AID PROCEDURES

May be harmful if swallowed, inhaled, or allowed to absorb through the skin. Wash contacted area with soap and water for 15 minutes. If inhaled remove to fresh air. Report exposure to the appropriate safety official. Consult a physician immediately.

SPECIAL HANDLING PRECAUTIONS

VENTILATION:	Appropriate ventilation to prevent the formation and dispersion of aerosols is
	required. When possible these materials should be handled in a designated
	biohazard work area that is segregated from common laboratory work areas.
EYE PROTECTION:	Required. Goggles or Safety glasses with side shields are recommended.
RESPIRATORY	Required. An approved respirator designed for use with toxic liquids is
PROTECTION:	necessary to prevent inhalation of aerosols.
PROTECTIVE GLOVES:	Required when handling any of these materials.
ADDITIONAL:	A full length laboratory gown with sleeves is required to prevent contact of any
	spilled liquids with clothing or exposed skin.

SPECIAL PRECAUTIONS

This material is for research and experimental application only. It is not intended for food, drug, household, agricultural, or cosmetic use. All materials should be handled only by technically qualified individuals experienced with working with potentially hazardous chemicals. The above information is correct to the best of our knowledge. The user should make independent decisions regarding completeness of the information, based on all sources available. EY Laboratories, Inc. shall not be held liable for any damage resulting from handling or contact with the above product.

EY LABORATORIES, INC. 107 North Amphlett Blvd. San Mateo, CA 94401

Tel:	650-342-3296
Fax:	650-342-2648
Orders:	1-800-821-0044
	(Outside CA only)